

# Fibertec™ 8000

The first fully automated fiber analysis solution following official reference methods



ANALYTICS BEYOND MEASURE

Fibertec™ 8000 is a fully automated system for determination of crude fibre, detergent fibre and related parameters according to standard reference 'crucible' methods such as Weende, van Soest etc.

### The lowest operator time of any fibre solution

Unattended measurement of up to six samples releases staff to do other things. On board heating and automatic dispensing of all reagents, antifoam and rinsing with water saves time and prevents the risk of human error.

### Official reference method results with unrivalled accuracy

No sample transfer or risk of error as sample residue remains in the crucible throughout the procedure. Discrete analysis of each sample according to official methods ensures reliable reference results, while consistent heating rate and time to boiling ensures repeatability.

### Safest fibre analysis solution available

Ensure safety with automatic functions such as dispensing of reagents to prevent contact with chemicals and fumes, heat reduction at boiling point to prevent spillage and addition of enzymes and anti-foam when required.

### Sample type

Raw materials and finished products in feed and agriculture

### Parameters

Crude Fibre (CF), Neutral Detergent Fibre (NDF), amylase treated Neutral Detergent Fibre (aNDF), Acid Detergent Fibre (ADF), and Acid Detergent Lignin (ADL)

### Technology

Fully automated 6 position single or sequential extraction including boiling, rinsing and filtration

### Official reference method

ISO 6865 & 92/89 EEC Crude Fibre in feed  
ISO 16472 aNDF in animal feed  
AOAC 2002.04 aNDF in feed  
ISO 13906 ADF and ADL in animal feed

# Specifications

| Performance data                                      |   |
|---|---|
| Sample size   | 0.5 - 3 g   |
| Measuring range                                       | 0.1% - 100%   |
| Capacity per batch                                    | Up to 6 samples simultaneously  |
| Capacity per day                                      | Up to 36 analyses (Crude Fibre method).<br>Up to 60 analyses using modified procedure |
| Repeatability   | ±1 % relative at 5% - 30% fibre level   |
| Reagent preheating time                               | 10 - 12 minutes   |
| Heating-up time from preheated temperature to boiling | 5 - 7 minutes   |

| Installation requirements:                                 |                   |                   |                         |        |  |
|--|-------------------|-------------------|-------------------------|--------|--|
| Equipment  | Power supply      | Power consumption | Dimensions w × d × h cm | Weight | Water supply   |
| Fibertec™ 8000 Hot Extraction Unit                         | 230 V<br>50/60 Hz | 2.000 W           | 73 × 39 × 64            | 67 kg  | *Tap water minimum 2 l/min (4-25°C, depending on water pressure) |
| FT 121 Fibertec™ Cold Extraction Unit with water aspirator | -                 | -                 | 56 × 38 × 28            | 14 kg  | Tap water 2 l/min  |

Note: Cold Extraction Unit must be placed in the fumehood with at least 0,5 m/sec airflow.

\* When Fibertec™ 8000 is in standby mode the tap water supply is closed down.

## System description:

Fibertec™ 8000, complete system, 230V, 50/60Hz comprising:

- Fibertec™ 8000, hot extraction unit
- FT 121 Fibertec™, cold extraction unit
- Standard accessory kit,
- Document kit

Fibertec™ 8000 system, 230V, 50/60Hz same as above but without cold extraction unit.

## Accessories:

Crucible stand for 6 crucibles, crucible holder, acid tank, alkali tank, NDS tank, ADS tank

## Optional accessories:

Crucibles, P0 (porosity 160 - 250 µm), set of 6  
Crucibles, P1 (porosity 100 - 160 µm), set of 6  
Crucibles, P2 standard (porosity 40 - 100 µm), set of 6  
Crucibles, P2 US (porosity 40 - 60 µm), set of 6  
Crucibles, P3 (porosity 16 - 40 µm), set of 6

000 «Диаэм»

Москва  
ул. Магаданская, д. 7, к. 3 ■ тел./факс: (495) 745-0508 ■ sales@dia-m.ru

www.dia-m.ru

С.-Петербург  
+7 (812) 372-6040  
spb@dia-m.ru

Новосибирск  
+7(383) 328-0048  
nsk@dia-m.ru

Воронеж  
+7 (473) 232-4412  
vrn@dia-m.ru

Йошкар-Ола  
+7 (927) 880-3676  
nba@dia-m.ru

Красноярск  
+7(923) 303-0152  
krsk@dia-m.ru

Казань  
+7(843) 210-2080  
kazan@dia-m.ru

Ростов-на-Дону  
+7 (863) 303-5500  
rnd@dia-m.ru

Екатеринбург  
+7 (912) 658-7606  
ekb@dia-m.ru

Кемерово  
+7 (923) 158-6753  
kemerovo@dia-m.ru

Армения  
+7 (094) 01-0173  
armenia@dia-m.ru

