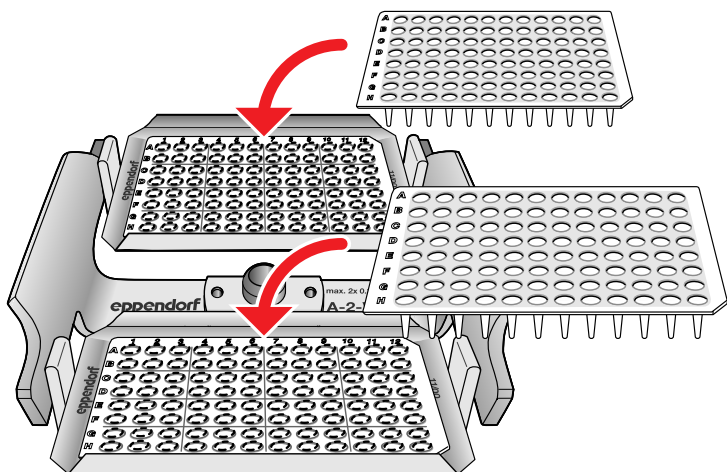
Mikrotiterplattenrotor A-2-VC / Microtiter plate rotor A-2-VC

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Unskirted PCR plates, PCR strips or single PCR tubes should be inserted into the Eppendorf work tray (order number: 0030 124.235) prior to centrifugation.

The work tray fits into the rectangular opening in the bottom of the bucket. There are guide slots which provide assistance with inserting the tray!

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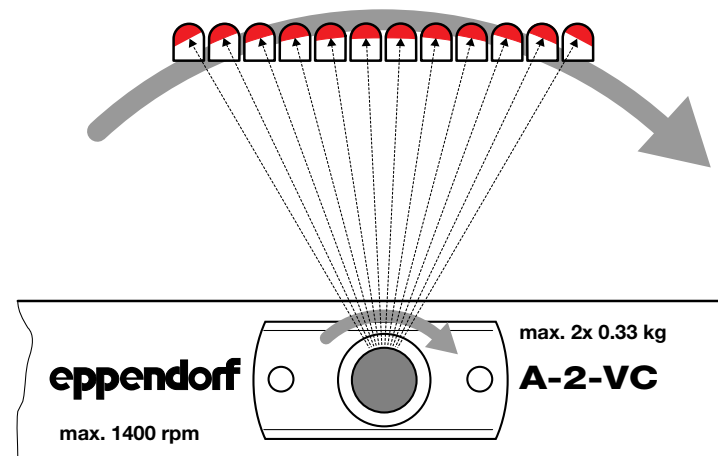
Please note: The liquid levels in the wells of the plate are perpendicular to the rotor axis during centrifugation. Due to the centrifugal force, the liquid levels in the wells at the edges of the plate are sloped.

 **This behavior can cause an over-flow if the liquid volume in the wells is too high!**

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Note:

When fully loaded (two microtiter plates per bucket), it may be possible to open lid of the Vacufuge before the rotor has come to a standstill.

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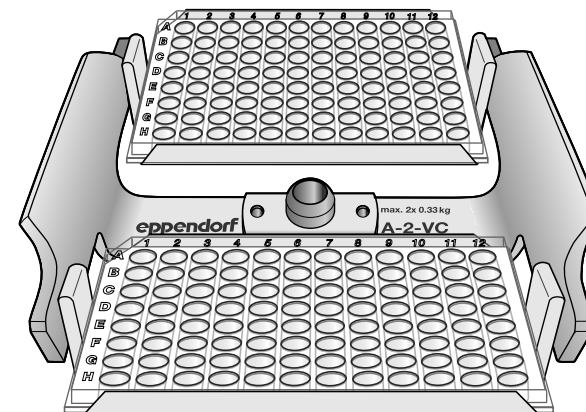
eppendorf

pendorf is a registered trademark.
90 045.963-01/0206 · Printed in Germany · Printed on chlorine-free bleached paper

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The following plates can be inserted directly into the microtiter plate buckets of the rotor:

Microtiter plates (max. two per bucket) or skirted PCR plates.

The plates must fit securely into the bucket. Avoid placing them partially outside of the bucket. Make sure that the plates are inserted into the bucket correctly. For safety reasons, please carry out a manual swing test with empty plates is recommended.

! Incorrectly positioned plates can cause damage!

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